



The Golden Gate™ Low Temperature Diamond ATR System



A variable temperature diamond ATR system for the FTIR measurement of solids and liquids at ambient to near liquid nitrogen temperatures.

The Golden Gate low temperature diamond ATR system is the first ATR accessory to provide high performance ATR measurements down to near liquid nitrogen temperature.

The system uses a thermally insulated copper and stainless steel dewar in conjunction with an integral heater to provide continuous temperature control across the working range. The high thermal conductivity of the diamond crystal provides rapid temperature stabilization, accurate temperature measurement, and avoids temperature gradients across the sample.

The diamond ATR crystal is high temperature metal bonded into a tungsten carbide support disk and the top plate is hard-anodized to make the accessory chemically resistant and capable of withstanding the pressures required for optimum crystal and sample contact. Pressure is applied to the sample using the quick lock and release bridge from the proven Golden Gate diamond ATR system. This clamping device allows rapid sample throughput and reproducible solid sampling. It also has a built-in torque limiter to control the loads applied to the diamond.

Key Sampling Benefits

- Continuous control of sample temperature from near liquid nitrogen temperatures to 80 °C.
- High thermal conductivity provides rapid cooling and temperature stabilization.
- Proven, strong clamping device, based on the Golden Gate Diamond ATR, allows rapid, reproducible sample throughput.
- Thermally insulated copper and stainless steel dewar allows for the use of liquid nitrogen, dry ice, or salt and water mixtures.
- Permanently loaded diamond ATR crystal is self-aligning, eliminating the need for optical re-alignment after cooling.

The Golden Gate low temperature diamond ATR is constructed in such a way that the crystal mounting is under a constant load. This ensures that the diamond is kept in constant optical alignment, negating the effects of thermal expansion and contraction. The upper body of the system is separated from the top-plate by a replaceable, thermally conducting spacer, avoiding the need for use of sealants that may contaminate the diamond surface. The system is easily and quickly taken apart for cleaning.

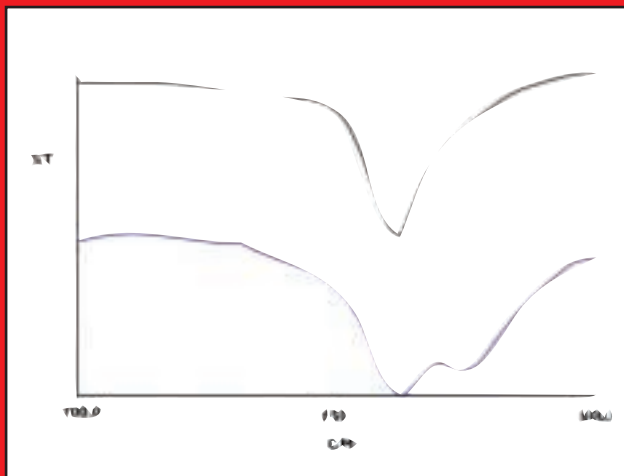


Figure 1: PTFE recorded at room temperature (black) and at -100 °C (blue) using a DTGS detector at 4 cm⁻¹ resolution.

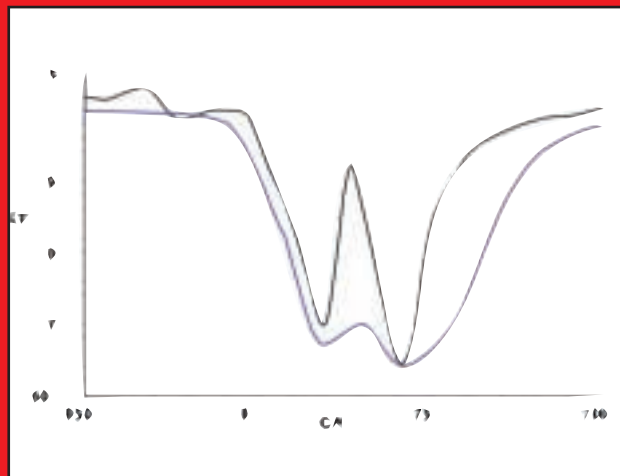


Figure 2: Carbon tetrachloride measured at room temperature (blue) and -143 °C (black) using a DTGS detector at 4 cm⁻¹ resolution.

Ordering Information

10592 Golden Gate Low Temperature Diamond ATR System, includes optics unit with choice of ZnSe or KRS -5 lenses, low temperature diamond top plate and temperature controller, sample clamping system with integrated torque limiter, Benchmark™ Instrument Base-plate and purge bellows.

Please specify instrument make and model, required voltage and country of destination.

10590 Golden Gate Low Temperature Diamond ATR Top Plate with temperature controller only, includes sample clamping system with integrated torque limiter.

Please specify instrument make and model, required voltage and country of destination.

Spares and Consumables

- 10593** Glass filled PTFE gaskets (pkt 5)
- 10594** High thermal transfer gaskets (pkt 5)
- 10595** Liquid sample injection device

Illustrations, descriptions and specifications - This Data Sheet was correct at the time of going to print. However, Specac's policy is one of continuous product development. The right is reserved to change descriptions and specifications at any time. For the latest details please contact your local Specac Representative. Golden Gate is a trademark of Specac Ltd.

smiths
A part of Smiths Group plc

Specac



SPECAC LIMITED,
River House, 97 Cray Ave, St. Mary Cray,
Orpington, Kent BR5 4HE, U.K.
Tel : +44 (0)1689 873134
Fax : +44 (0)1689 878527
www.specac.com
Registered No. 1008689 England

SPECAC INC.,
410 Creekstone Ridge,
Woodstock,
GA 30188, USA
Toll Free : +1 800 447 2558
Fax : +1 770 319 2488
Specac is part of Smiths Group plc